

TenuPol-5



Automatic electrolytic thinning of specimens for transmission electron microscopy



Unique advantages and technologies

Scan function

Determination of correct polishing voltage through built-in scan function

Built-in database

With 18 Struers and up to 200 user methods

Automatic shut-off

Thinning process stopped automatically through infra-red light when hole appears

Pre-thinning

Large samples can be pre-thinned electrolytically to avoid mechanical deformation


Simultaneous thinning

Thinning from both sides to avoid damaged structures.


With TenuPol-5 a perforated specimen for transmission electron microscopy can be made from a sample of 3 or 2.3 mm dia. in just a few minutes.

The specimen is polished from both sides simultaneously, thus providing a structure with a minimum of deformation. When the perforation appears, the polishing can automatically be stopped by the infrared detector system, and the specimen is ready for the TEM examination. No special operator training is required.


Also electrolytical pre-thinning or blanking can easily be carried out.

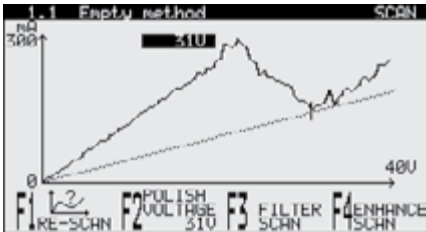

Determination of polishing voltage through built-in scan function
TenuPol-5 is equipped with a unique scanning function. After placing a specimen in the holder, a pre-defined voltage range is scanned to determine the current density curve. This curve is used to define the correct polishing voltage for the thinning process. Instead of time consuming trial and error testing of parameters it is now easy to find the correct settings after a single scan of a new sample material.

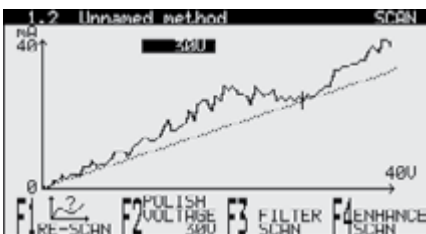

As not all materials display such an easy to analyse current density curve as shown before, both filter and enhance functions are included with the TenuPol-5.

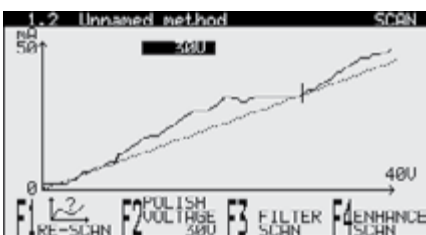

After filtering the current density curve might look like that, and the correct polishing value is easy to define.



After an additional enhancement the definition might be even easier.

STRUERS METHODS			
1	Material	Polishing Voltage	Filter/Enhance
1	Stainl. steel	0300V	0 *
2	Low C-steel	0100V	0 *
3	Low C-steel	0200V	0 *
4	Apax 45HR	0100V	0 *
5	Apax 45HR	0200V	0 *
6	Apax 45HR	0300V	0 *
7	Titanium	0100V	0 *
8	Titanium	0200V	0 *

Built-in database with 18 Struers and up to 200 user methods

10 methods for the final thinning and 8 methods for pre-thinning of various materials are included with TenuPol-5. A whole range of materials can thus be prepared immediately, without any lengthy and time consuming trials. These methods can also be used as starting point for the development of methods for other materials. 200 user-defined methods can be saved in 20 groups in the database of TenuPol-5.

1.1 Stainl. steel 0300V EDIT METHOD	
Electrolyte :	08
Voltage :	50.00
Temperature recomb. :	+15°C (-5.0°C)
Polishing time :	No limit
Light stop value :	Auto
Graphic time scale :	Auto
Flow mode :	Single flow
Pump flow rate :	12

Automatic shut-off

The very small perforations required for the following examination in the TEM are very difficult to control visually.

TenuPol-5 is therefore equipped with an IR-detector system. The infrared

light from the transmitter passes through the jet nozzle to the sample. As soon as a hole appears in the centre of the specimen, the light hits the photocell on the other side and the process is shut off in accordance with the pre-set light stop value. As the IR-detector system is insensitive to white light, errors are eliminated.

When pre-thinning or "punching" specimens, the automatic shut-off system using infrared light cannot be used. Then a preparation time is set instead.

Pre-thinning

To avoid deformations of the finished sample the basic material with a diameter of 12-21 mm is pre-thinned to a thickness of less than 0.5 mm in an area of up to 10 mm dia. This procedure is carried out by means of a special specimen holder.

Electrolytic "punching"

Mechanical punching to produce samples small enough for electrolytical thinning is bound to introduce deformations in the sample material. With TenuPol-5 several 3 mm (or 2.3 mm) specimens at a time can be "punched" electrolytically from the foil material, without introducing any deformation at all.


Chemical thinning

TenuPol-5 is made of corrosion-resistant materials. This allows for chemical polishing with most chemicals. TenuPol-5 can be used with perchloric acid and other highly corrosive electrolytes.

Design

TenuPol-5 consists of two separate units, the control unit and the polishing or thinning unit.

The polishing unit with the electrolyte pump, electronic thermometer and exchangeable specimen holder

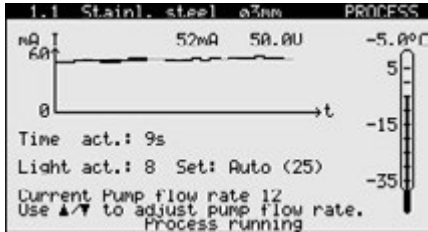

The control unit

Both the power supply and the programming and monitoring functions are incorporated in the control unit. The unit is operated on the touch pad and on the backlit graphic display all parameters are presented. Upon pressing the Start button, the display changes, and the ongoing process can be monitored. Parameters like current, electrolyte temperature and elapsed thinning time are displayed. Both the display of parameters before and the display of actual values after process start provide an exceptional amount of immediately usable information. Deviations from one process to the next can be detected straightaway, and necessary measures can be taken. Required changing of electrolyte can be detected before changes of the polishing result will be noticed.

The polishing/thinning unit

The design as a separate unit makes it possible to place the polishing/thin-

ning unit away from the control unit. If required the polishing unit can be set up in a fume cabinet while the control unit is set up outside. TenuPol-5 comprises 2 electrolyte reservoirs, an insulated one for use with the built-in cooling coil and an uninsulated one for cooling the electrolyte, e.g. by placing the reservoir in an ice bath.

Specimen holders

The TenuPol-5 standard equipment includes a specimen holder for 3 mm specimens. A 2.3 mm specimen holder and a sample holder with an opening of 10 mm dia. are also available. The holders are made of reinforced Teflon and consist of two main parts. The holders are easily opened and closed, allowing simple mounting and fast cleaning. One of the two parts carries the platinum conductor. When the closed holder is mounted in the polishing cell, electrical connection to the polishing

circuit is automatically established. No extra plug or cable is necessary. The other part has a sliding diaphragm which allows the mounting of samples with any thickness less than 0.5 mm. The 3 mm and 2.3 mm specimen holders are normally used for final thinning (perforation), and the holder with a 10 mm diaphragm for pre-thinning and electrolytic "punching".

Jet holders

The standard set of jet holders has a 1 mm jet bore for use with the 3 and 2.3 mm specimen holders.

A set of jet holders with a 2.5 mm jet bore for use with the holder with a 10 mm opening is also available.



A specimen holder for 3 mm samples for final thinning, with 10 mm dia. hole for pre-thinning or electrolytic punching and a set of jet holders. The platinum connector and the sliding specimen fixture can be seen.

Technical data

Control unit

Mains voltage:	Single-phase, 100-120/220-240V, 50-60 Hz.
Power consumption:	220-240V / 1.50A 110-120V / 3.15A
Output:	0-100V / 0-2A

Dimensions

Width:	385 mm
Depth:	350 mm
Height:	160 mm
Weight:	17 kg

Polishing unit

Sizes of original specimen:	12-21 mm dia., max. 1 mm thick 3 mm dia., max. 0.5 mm thick 2.3 mm dia., max. 0.5 mm thick
-----------------------------	--

Dimensions

Width:	270 mm
Depth:	180 mm
Height:	276 mm
Weight:	3.8 kg

Specifications

	Code
TenuPol-5 control unit, with electronic thermometer and adapter for the connection of (04086001)	05396133
TenuPol polishing unit, with specimen holder for 3 mm dia. specimens (04086901), sets of jets (04086904), pump, cooling coil, insulated PVC container, non-insulated PVC-container and built-in photo cell	04086001
Specimen holder for 3 mm dia. specimens	04086901
Specimen holder for 2.3 mm dia. specimens	04086902
Specimen holder, 10 mm diaphragm for pre-thinning	04086903
Set of jets, 1 mm bore	04086904
Set of jets, 2.5 mm bore for use with (04086903)	04086905
Tape Kit for electrolytic "punching"	40900041

Struers' equipment is in conformity with the provisions of the applicable International Directives and their appurtenant Standards. (Please contact your local supplier for details).

Struers' products are subject to constant product development. Therefore, we reserve ourselves the right to introduce changes in our products without notice.



Struers A/S

Pederstrupvej 84
DK-2750 Ballerup, Denmark
Phone +45 44 600 800
Fax +45 44 600 801
struers@struers.dk
www.struers.com

AUSTRALIA & NEW ZEALAND

Struers Australia
27 Mayneview Street
Milton QLD 4064
Australia
Phone +61 7 3512 9600
Fax +61 7 3369 8200
info.au@struers.dk

BELGIQUE (Wallonie)

Struers S.A.S.
370, rue du Marché Rollay
F- 94507 Champigny
sur Marne Cedex
Téléphone +33 1 5509 1430
Télécopie +33 1 5509 1449
struers@struers.fr

BELGIUM (Flanders)

Struers GmbH Nederland
Elektraweg 5
3144 CB Maassluis
Telefoon +31 (10) 599 7209
Fax +31 (10) 5997201
netherlands@struers.de

CANADA

Struers Ltd.
7275 West Credit Avenue
Mississauga, Ontario L5N 5M9
Phone +1 905-814-8855
Fax +1 905-814-1440
info@struers.com

CHINA

Struers Ltd.
No. 1696 Zhang Heng Road
Zhang Jiang Hi-Tech Park
Shanghai 201203, P.R. China
Phone +86 (21) 6035 3900
Fax +86 (21) 6035 3999
struers@struers.cn

CZECH REPUBLIC

Struers GmbH
Organizační složka
Havlíčková 361
CZ-252 63 Roztoky u Prahy
Phone +420 233 312 625
Fax +420 233 312 640
czechrepublic@struers.de

DEUTSCHLAND

Struers GmbH
Carl-Friedrich-Benz-Straße 5
D-47877 Willlich
Telefon +49 (0) 2154 486-0
Fax +49 (0) 2154 486-222
verkauf@struers.de

FRANCE

Struers S.A.S.
370, rue du Marché Rollay
F-94507 Champigny
sur Marne Cedex
Téléphone +33 1 5509 1430
Télécopie +33 1 5509 1449
struers@struers.fr

HUNGARY

Struers GmbH
Magyarországi Fióktelep
Tatai ut 53
2821 Gyermely
Phone +36 (34) 880546
Fax +36 (34) 880547
hungary@struers.de

IRELAND

Struers Ltd.
Unit 11 Evolution@ AMP
Whittle Way, Catcliffe
Rotherham S60 5BL
Tel. +44 0845 604 6664
Fax +44 0845 604 6651
info@struers.co.uk

ITALY

Struers Italia
Via Monte Grappa 80/4
20020 Arese (MI)
Tel. +39-02/38236281
Fax +39-02/38236274
struers.it@struers.it

JAPAN

Marumoto Struers K.K.
Takara 3rd Building
18-6, Higashi Ueno 1-chome
Taito-ku, Tokyo 110-0015
Phone +81 3 5688 2914
Fax +81 3 5688 2927
struers@struers.co.jp

NETHERLANDS

Struers GmbH Nederland
Elektraweg 5
3144 CB Maassluis
Telefoon +31 (10) 599 7209
Fax +31 (10) 5997201
netherlands@struers.de

ÖSTERREICH

Struers GmbH
Zweigniederlassung Österreich
Betriebsgebiet Puch Nord 8
5412 Puch
Telefon +43 6245 70567
Fax +43 6245 70567-78
austria@struers.de

POLAND

Struers Sp. z o.o.
Oddział w Polsce
ul. Jasnogórska 44
31-358 Kraków
Phone +48 12 661 20 60
Fax +48 12 626 01 46
poland@struers.de

ROMANIA

Struers GmbH
Sucursala Sibiu
Str.Scoala de Inot, nr. 18
RO-550005 Sibiu
Phone +40 269 244 558
Fax +40 269 244 559
romania@struers.de

SCHWEIZ

Struers GmbH
Zweigniederlassung Schweiz
Weissenbrunnstraße 41
CH-8903 Birmensdorf
Telefon +41 44 777 63 07
Fax +41 44 777 63 09
switzerland@struers.de

SINGAPORE

Struers Singapore
627A Aljunied Road,
#07-08 BizTech Centre
Singapore 389842
Phone +65 6299 2268
Fax +65 6299 2661
struers.sg@struers.dk

SUOMI

Struers Suomi
Kalevankatu 43
00180 Helsinki
Puhelin +358 (0)207 919 430
Faksi +358 (0)207 919 431
finland@struers.fi

SWEDEN

Struers Sverige
Ekbacksvägen 22
168 69 Bromma
Telefon +46 (0)8 447 53 90
Telefax +46 (0)8 447 53 99
info@struers.se

UNITED KINGDOM

Struers Ltd.
Unit 11 Evolution @ AMP
Whittle Way, Catcliffe
Rotherham S60 5BL
Tel. +44 0845 604 6664
Fax +44 0845 604 6651
info@struers.co.uk

USA

Struers Inc.
24766 Detroit Road
Westlake, OH 44145-1598
Phone +1 440 871 0071
Fax +1 440 871 8188
info@struers.com